
PRS-2000 Compact Protein Review Station

FOR X-RAY CRYSTALLOGRAPHY

STREAMLINING THE TASK OF IDENTIFYING PROTEIN CRYSTALS — WITH ASTONISHING NEW BREAKTHROUGHS IN COST AND CONVENIENCE

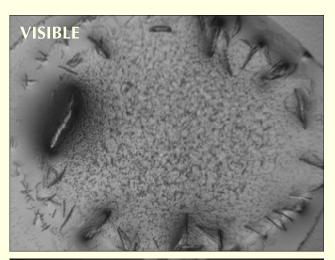



Korima's Ultraviolet Technology Revolutionizes Protein Crystal Recognition

- Fail-safe protein crystal identification Every time—in every sample
 - Extremely small footprint
 - Easy to operate
 - Quick setup
 - New low price

Traditionally, X-ray crystallographers have faced a truly daunting challenge—that of scanning hundreds of potentially successful images for protein crystal growth.

Even though microplates of buffer solutions could be teeming with critical protein crystals, target proteins often could go undetected among strong salt growth and other debris.

Experience, of course, would help a crystallographer discern target proteins. But with so many decoys appearing in countless shapes and sizes, it often took a measure of sheer luck to pick out the crystals that really counted.

ENTER THE PRS-2000

The PRS-2000 Protein Review Station instantly revolutionizes the entire protein recognition process.

Under Korima's exclusive, patentpending split fluorescent beam, target proteins—and *only* target proteins—respond selectively to fluorescent excitation.

With the PRS-2000, you will instantly recognize protein crystals. All of them—in every solution. Eliminating guesswork. Saving time. Slashing costs.





PRS-2000

Features and **Specifications**

Breakthrough Technology that Fits in Anywhere!

The PRS-2000 makes a major difference in any pharmaceutical or research lab—with the smallest footprint ever!

Fitting easily on a standard lab bench, the complete system comprises all necessary components, including microscope, dual cameras and monitors, lights, and power supplies.

- LCD Monitors. 15- and 19inch LCD monitors display samples simultaneously under ultra-violet and visible light.
- Microscope and Objectives.
 Turret-mounted 5x, 10x, and 20x UV/Visible objectives enable complete cell coverage or close inspection of individual proteins.
- PC and Software. A dedicated Pentium IV/Windows XP Professional PC and proprietary software stand at the heart of the system.



The PRS-2000 Workstation fits easily on a lab bench.

- Lights and Cameras. Visible and shutter-controlled DUV cameras capture the effects of highcontrast visible and ultra-violet lighting.
- X-Y Platform. The system's standard X-Y platform allows precise, systematic examination of samples in a 96-well Limbro plate or any standard microplate.

An optional, motorized X-Y table, which allows automated archiving, is also available.



119 E. Star of India Lane Carson, CA 90746 www.korimainc.com Voice: (310) 532-2222 Fax: (310) 527-0470

email: info@korimainc.com